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型式 LA-150CE
LA-150UE-B
LA-150UE-C

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16 WARRANTY AND AFTER-SALE SERVICES

• A warranty period for this product is one year counting from the date of purchase. During this period, in spite of using the product in normal operating conditions, if you encounter any trouble, we will repair it free from charge.

I. Troubles or damages attributable to wrong operation, re-engineering, or improper repair.
II. Troubles or damages attributable to drop, or transportation of the product after purchase.
III. Troubles or damages attributable to natural disasters such as fire, earthquake, flood disaster, lightning, and environmental pollution, abnormal voltage, and other external factors.

IV. When the warranty card is not presented. (use the warranty card separately provided.)
V. When the date of purchase, customer's name, and dealer's name are not entered in the warranty card, or when the words and purchase are rewritten.

VI. Replacement of consumable parts. (halogen lamp, lamp socket, etc)

※ This warranty is applicable only domestically. For the troubles resulting from overseas use, consult the dealer where you bought the apparatus or our office.

◎ When you ask for a repair, please let us know the following :

- Model number :
 - Manufacturer number :
 - Details of trouble
 - Date of purchase
 - Your name, company name, and section
 - Address, phone number / fax number
- the apparatus.
number sticker attached to
inscribed on the manufacture

HAYASHI - REPIC CO., LTD

Division-1 Office : 1-28-3, Kitaohtsuka, Toshima-ku, Tokyo
Phone : 03 - 3918 - 5623 FAX : 03 - 3918 - 5683
Sendai Sales Office : 2-8-8, Uesugi, Aoba-ku, Sendai-shi, Miyagi-ken
Phone : 022 - 221 - 0471 FAX : 022 - 221 - 0472
Osaka Sales Office : 9 - 12, Hirotsuba-cho, Suita-shi, Osaka-fu
Phone : 06 - 6369 - 5023 FAX : 06 - 6369 - 5022

13. MAINTENANCE PARTS

- Consumable parts such as halogen lamp, lamp socket, fuse, are in stock as spare parts. When you want to procure the maintenance parts, contact our dealer where you purchased this product or our office.
- Halogen lamp : • JCR15V150W 5H/5 • JCR12V100W10H • JCR12V100W10H
- Lamp socket : • HS-310-100-S (Life: 3,000 hours)
- Fuse : • 250V T4A

14. TROUBLE SHOOTING (Check again before asking for a repair)

Trouble	Check	Remedy
• The light source dose not work even if the power is turned on. (Lamp unilluminated. Fan not running.)	• Whether or not the Power plug is connected to the plug socket. • Whether or not the fuse is gone. • Whether or not the overheat protective circuit is activated.	• Connect the Power plug to the plug socket. • Replace the fuse. • Eliminate a cause for high temperature, cool down the light source, and then, turn on the power switch again.
• Even if the power is turned on, the lamp is not illuminated; (Fan running, but lamp unilluminated.)	• Whether or not the lamp is burnt out. • Whether or not the lamp socket is worn out. • Whether or not the light quantity control knob is set to 0.	• Replace the lamp. • Replace the lamp socket. (Life : 3,000 hours) • Turn the light quantity control a light quantity.
• A light quantity dose not be changed by turning the light quantity control knob.	• [MANUAL / REMOTE] selector switch is set properly.	• Set the switch to [MANUAL]
• Light Intensity Remote function doesn't work.	• Whether or not the [MANUAL / REMOTE] selector switch is set properly. • Whether or not the light intensity remote control voltage is impressed. • Whether or not the noise ripple touches the pin of the light intensity remote control voltage. • Whether or not there is plenty of current offered to the unit which provide voltage. (prepared by customer)	• Connect to the unit which provide voltage 0-5VDC • Take measures with noise ripple to the unit and connect cable which prepared by customer. • Make sure the current is equal to which mentioned in [EXTERNAL CONTROL].

If the product has troubles, unplugging the power supply-cord immediately to disconnect the power. Please be sure to contact the manufacturer for repair.


15. CHANGING THE FUSE

When the fuse exchanged, unplugging the power supply-cord to disconnect the power.


• How to Replace the Fuse

- ① Lightly pushing a fuse holder's cap with a flat screwdriver, turn in to the left to unlock the cap.
- ② Pull out the cap from the fuse holder. Remove a snapped fuse from the cap.
- ③ Set a replacement fuse into the cap. Put back the cap into a fuse holder.
- ④ Lightly pushing the cap with a flat screwdriver, turn in to the right to lock the cap.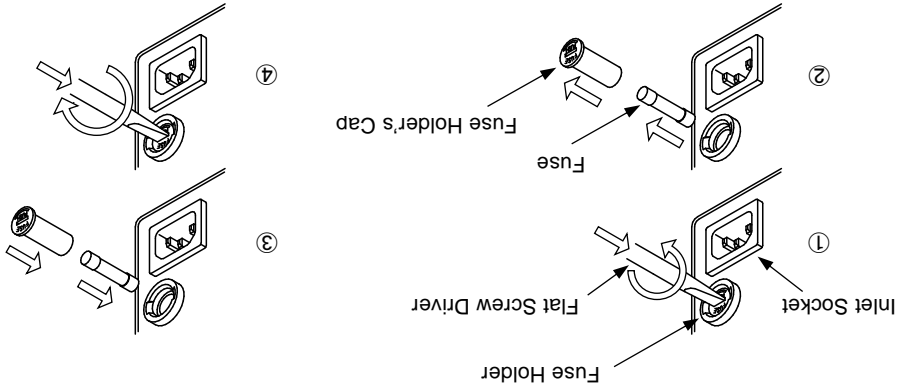


11. EXTERNAL DIMENTION DRAWING

LA-150UE-B, LA-150UE-C Dimensions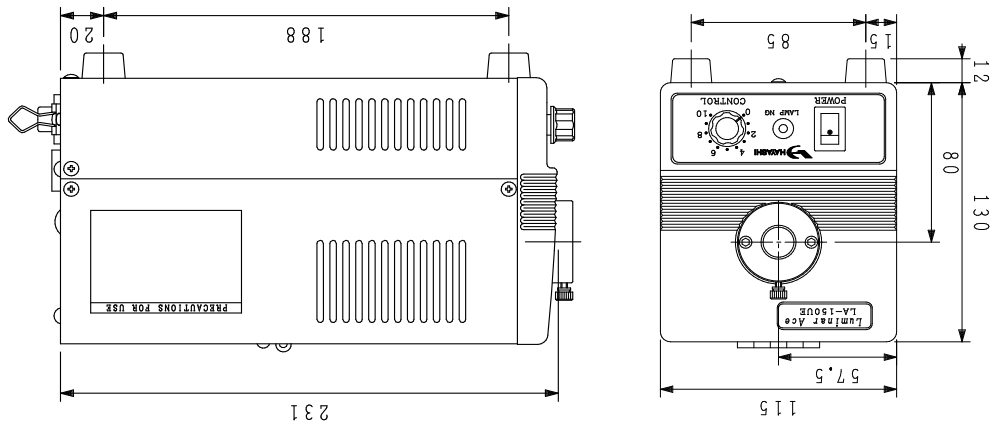

LA-150CE Dimensions

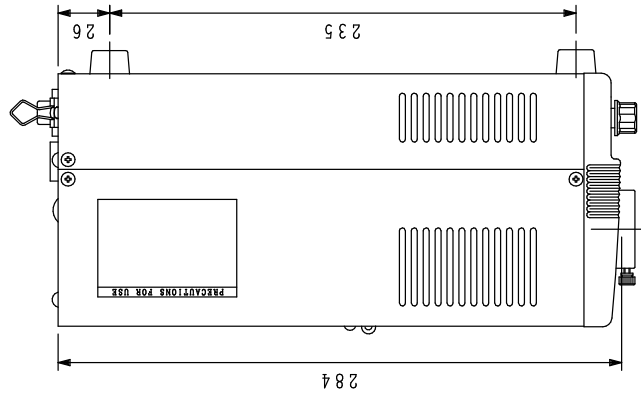12. LIGHT GUIDE

Precautions for Use

1) The light guide is a bundle of fine glass fibers. Note that if it is pulled hard or bent beyond allowable limits, the fibers will snap.

2) When the light guide is badly contaminated, wipe it quickly with alcohol.

Never use solvent such as thinner, benzine, acetone, and so on. They may dissolve the product.

ATTENTION







When you use a light guide, be sure to use the item specified by the manufacturer. For further information, please inquire to the manufacturer.

10. SPECIFICATION

(The specifications of the product are subject to change without prior notice due to its improvement.)

Model number	LA-150CE	LA-150UE-B	LA-150UE-C
*1 Input voltage	200 to 240VAC±10%	100 to 120VAC±10%	200 to 240VAC±10%
Rated frequency	50 / 60Hz		
Power consumption	180W (At 150W lamp maximum rated load)		
Operating	15V / 12V lamp available (With output voltage selector)		
Halogen lamp types	JCR15V150W5H/5 / JCR15V150WB Color temperature : 3,100K – 3,200K *2 Average life : 700hours < 15V halogen lamp > JCR12V100W10H / JCR12V100WB Color temperature : 3,000K – 3,100K *2 Average life : 1,000hours JCR12V50W20H Color temperature : 3,000K *2 Average life : 2,000hours		
Lighting system	DC lighting system		
Light quantity control	Output voltage variable system (With output selector) Variable voltage range : 3 to 12V / 14VDC Light quantity variable range : 2 to 100%		
Protective functions	Overcurrent protective circuit : Droop type (Automatic reset system) Overheat protective circuit : Output shut-off type (Reset by turning on the power again) Forced cooling by fan		
Cooling method	Optical filter, Heat absorbing filter incorporated		
Operating fuse	250VAC, T4A glass tube fuse (φ5.2 x 20mm)		
External control	① Lamp ON / OFF function		
Function	② Light quantity remote control function		
Other Function	Lamp burnout detect control (With monitor LED)		
Power supply-cord	For Japan: PSE-listed power supply-cord. For countries excluding Japan where the power supply is 100 to 120VAC: UL-listed power supply-cord. For countries where the power supply is 220 to 240VAC: Use the power supply-cord specified in the countries.		
Operating conditions	Temperature: 0°C to 40°C Humidity: 20% to 80%RH (no condensation) Pollution degree: Degree 2 Altitude: 2,000m max. Overvoltage category: Category II Short-term temporary overvoltage: 1,440V (up to 5 seconds) Long-term temporary overvoltage: 490V (longer than 5 seconds) Indoor use only		
Storage conditions	Temperature: -20°C to 60°C, Humidity: 90%RH max. (no condensation) UL, cUL, CE UL, cUL, PSE UL, cUL		
Safety	UL, cUL, CE		
Dimensions *3	W115×H130×D284 (mm)		
Weight	3.5kg		
	W115×H130×D231 (mm)		
	2.6kg		

*1 Input voltage: 150CE, 150UE-C "200-240Vac only", 150UE-B "100-120vac only",
 *2 Average life : The average lamp life is only information for customer, not a guarantee value.
 *3 Dimensions : Projections are excluded.
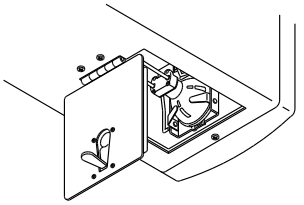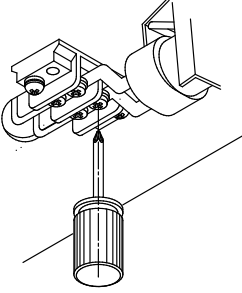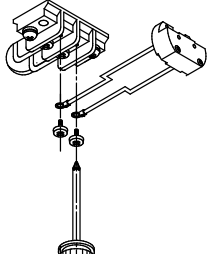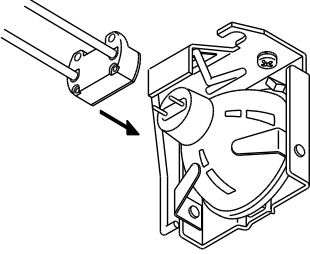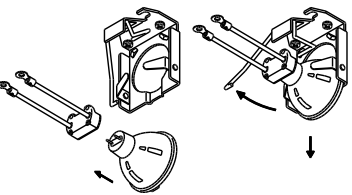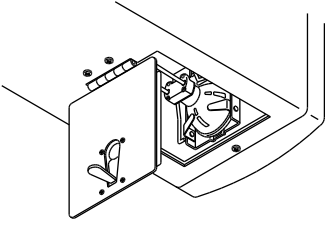
<p>① Release the cover stopper attached to the lamp house opening / closing cover on the top of this light source to fully open the cover.</p> 	<p>② Loosen the terminals board fixing screw and remove it.</p> 	<p>④ Fix the terminal of lamp socket to the terminals board with a screw.</p>  <p>[ATTENTION] Be sure to tighten the screw, if the screw doesn't be tightened, it may cause a failure.</p>	<p>⑤ Fix the lamp to the lamp holder</p> <p>Align the lamp terminals with the lamp socket and fix from in slowly.</p> 	<p>③ Operating the lamp removing lever of the lamp holder, push up the lamp. Hold the reflecting mirror of the lamp with fingers and draw out the lamp from the lamp holder.</p> 	<p>⑥ Close the cover slowly and fix it firmly with the cover stopper.</p> 
---	---	--	--	--	--

9. SHUTTER FUNCTION (OPTION)

- 9 - 1 Drive System
- The shutter used is of rotary solenoid drive system.

- 9 - 2 Driving Method

- Apply a voltage to between the no.1(+) and no.2(-) pins of an accessory shutter plug, the shutter will be activated.

The customer is requested to prepare the shutter drive power source. Use the power sources of the following capacities :

12VDC type : 12VDC, 1.0A or more
24VDC type : 24VDC, 0.5A or more


ATTENTION

Shutter OPEN / CLOSE

Drive Voltage of Solenoid	Type	SO type	SC type
0V	12 / 24V	close	open
		open	close

9 - 3 Shutter Receptacle

Shutter receptacle
• Receptacle type: HR10-7R-4S
• Adaptive Plug type: HR10-7P-4P (HIROSE)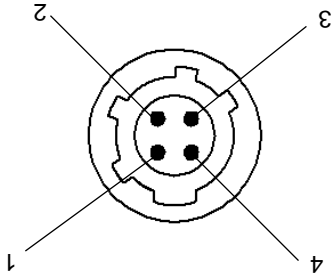

1 : Shutter drive supply voltage Input (+)
2 : Shutter drive supply voltage Input (-)
3,4 : Empty

8. Replacement of MAINTENANCE PARTS.









Be sure to use only specified lamp and lamp socket by manufacturer. Using of non-specified ones will lead to a trouble.

When the lamp and lamp socket exchanged, unplug the power supply-cord to disconnect the power.

Immediately after the lamp is turned off, it is reflecting mirror or lamp socket are still hot. To avoid burning your hand, wait until they are cooled down.

8 - 1 Lamp
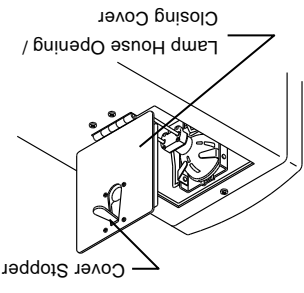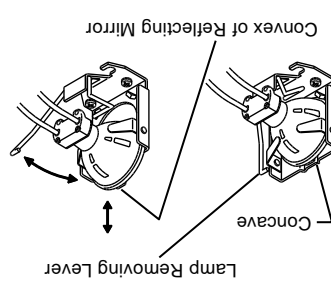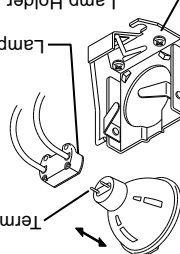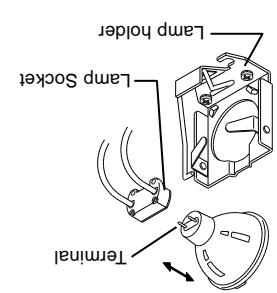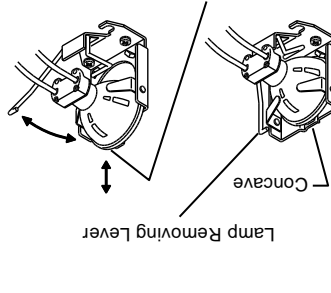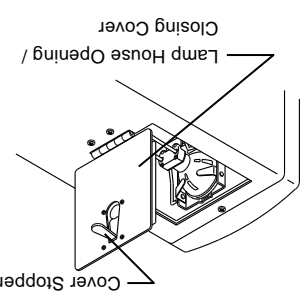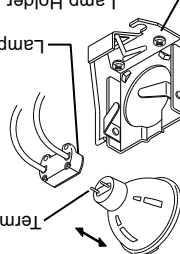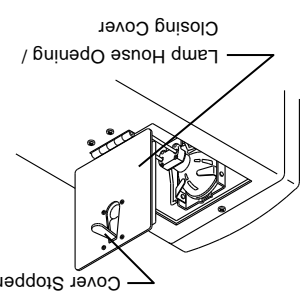
<p>① Release the cover stopper attached to the lamp house opening / closing cover on the top of this light source to fully open the cover.</p>  <p>Labels: Closing Cover, Lamp House Opening / Closing Cover, Cover Stopper.</p>	<p>② Operating the lamp removing lever of the lamp holder, push up the lamp. Hold the reflecting mirror of the lamp with fingers and draw out the lamp from the lamp holder.</p>  <p>Labels: Convex of Reflecting Mirror, Concave, Lamp Removing Lever.</p>	<p>③ Remove the lamp from the lamp socket.</p>  <p>Labels: Lamp Holder, Lamp socket, Terminal.</p> <p>INSPECTION</p> <p>When replacing the lamp, check the lamp socket for damages. When the lamp socket has been used for along period of time, the electrodes will be oxidized or corroded due to heat of the lamp or depending on the environment, and consequently, the service life of the lamp will be shortened. It is recommended to replace the lamp socket at least once a year.</p>	<p>④ Align the lamp terminals with the lamp socket and fit them in slowly.</p>  <p>Labels: Lamp holder, Lamp Socket, Terminal.</p> <p>ATTENTION</p> <p>Do not touch with the bare hand to the lamp bulb. If a lamp bulb is stained with fingerprints, etc., its service life may be reduced considerably or the lamp may burst.</p>
<p>⑤ Align the convex of the reflecting mirror of the new lamp with the concave of the lamp holder and push in the lamp into the lamp holder.</p>  <p>Labels: Convex of Reflecting Mirror, Concave, Lamp Removing Lever.</p>	<p>⑥ Close the cover slowly and fix it firmly with the cover stopper.</p>  <p>Labels: Closing Cover, Lamp House Opening / Closing Cover, Cover Stopper.</p> <p>ATTENTION</p> <p>Upon completion of lamp replacement work, turn on the power after confirm the opening / closing cover firmly closed.</p>	<p>⑦ Remove the lamp from the lamp socket.</p>  <p>Labels: Lamp Holder, Lamp socket, Terminal.</p> <p>INSPECTION</p> <p>When replacing the lamp, check the lamp socket for damages. When the lamp socket has been used for along period of time, the electrodes will be oxidized or corroded due to heat of the lamp or depending on the environment, and consequently, the service life of the lamp will be shortened. It is recommended to replace the lamp socket at least once a year.</p>	<p>⑧ Close the cover slowly and fix it firmly with the cover stopper.</p>  <p>Labels: Closing Cover, Lamp House Opening / Closing Cover, Cover Stopper.</p> <p>ATTENTION</p> <p>Upon completion of lamp replacement work, turn on the power after confirm the opening / closing cover firmly closed.</p>

7. EXTERNAL CONTROL


7 - 1 Remote Control Receptacle

Pin No.	I/O	Function	Internal Working condition	External Control System Connection
1	I	Input Lamp ON / OFF Signal (HOT)	12VDC pull up	Connect the switch (The allowed current should be more than 1.5mA) in lamp OFF for low level.
8	I	Input Lamp ON / OFF Signal (COLD)	GND level of second circuit	
2	I	Input intensity remote control voltage (HOT)	Allow voltage : 0 - 5VDC	Power supply for external control must meet the condition : low impedance output ; Voltage : less than 5 ; Current : more than 10mA ; NO noise ripple when connected.*1
9	I	Input intensity remote control voltage (COLD)	GND level of second circuit	
3	O	Output Lamp change signal (open collector)	Photo coupler output	Recommended use to control the circuit worked under the condition : less than DC 7mA, less than 24VDC*2
10	O	Output Lamp change signal (open emitter)		
4	I	11	2	empty
7	I	14	1	empty

- *1 It is used a shielded wire to prevent influence of outside noise, please shorten it as much as possible.(Recommended less than 2m)
 *2 Even though the lamp burning out signal doesn't output when the voltage reduce to less than 2.5VDC, it is not abnormal.


7 - 2 Receptacle

- Receptacle type: 57-40140
- Adaptive Plug type: 57-30140 (DDK Ltd.)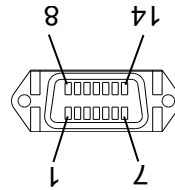

- 1-8: Halogen lamp ON / OFF
- 2-9: Light quantity Remote control voltage input (0 to 5VDC)
- 3: Lamp burnout signal output (open collector)
- 10: Lamp burnout signal output (open emitter)


7 - 3 Control Voltage Output Voltage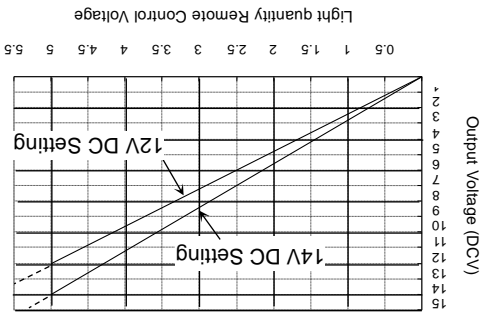

[ATTENTION] A remote control voltage should not be up to +5VDC at maximum.

If this limit is exceeded, an output voltage will go beyond the rating, leading to lamp burnout or power source failure.

7 - 4 External Control System Connection Diagram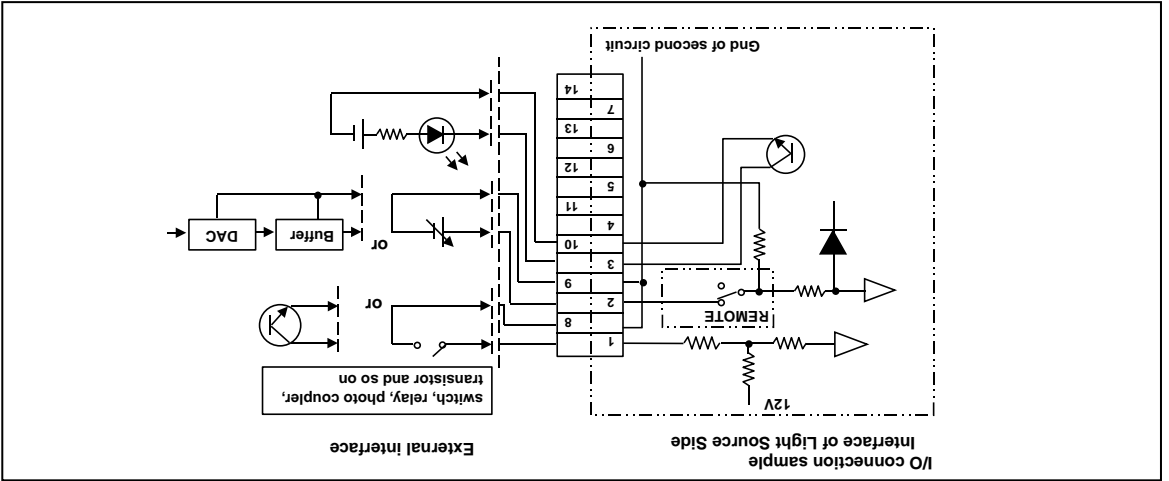

6. HOW TO USE

ATTENTION

If the product is used by the method which the manufacturer does not specify, the protection performance which product has is spoiled.

When you connect the remote control receptacle ⑪ or operate the 「MANUAL / REMOTE」 selector switch ⑩ or output voltage selector switch ⑧, be sure to turn off the power switch ① first in order to protect the power circuit in this product.

Be fully carefully not to confuse the set value of the output voltage selector switch with the rated voltage of the halogen lamp, because the following abnormalities will occur.

Output Voltage Selector Switch	Rated Voltage	Abnormality
12V	15V	• A light quantity (illumination) at maximum light quantity control will drop.
14V	12V	• The filament of the halogen lamp snaps the moment the power is turned on or the lamp life will be shortened considerably.

- 1) According to the lamp you want to use, set the output voltage selector switch ⑧ to 14V or 12V.
 Lamp used • 14VJ : 150W type
 • 12VJ : 100W / 50W type

- 2) According to the operation mode you want to choose, MANUAL or REMOTE mode.
 • 「MANUAL」 : Manual light quantity control (Use control knob)
 • 「REMOTE」 : Remote light quantity control (Use External control)

When using the apparatus at REMOTE, connect respective external control systems to an accessory remote control connector plug, and then, to the remote control receptacle ⑪. Refer to 「6.External」.

- 3) Insert the light guide in to the light guide socket ④ of the light source.

- 4) Adjust the light intensity to the minimum.

- 5) Connect the AC inlet ⑬ of the light source by power supply cord (100 – 240VAC line), turn on the power switch ①, and confirm that the apparatus is in a normal state.
 a) Make sure that the lamp burnout warning LED (red) ② is unilluminated.
 b) Make sure that the cooling fan in the exhaust port ⑦ is running.

- 6) After the lamp is turned on, set to your desired brightness.

< When manually controlling the light quantity >

Turn the light quantity control knob slowly (The light is getting brightness

by turning in the clockwise direction.)

> When remotely controlling the light quantity >

Apply a voltage, which can reach your desired light brightness to a light quantity remote control voltage input terminal.

(A remote control voltage should be applied within a range of 0 to 5 V DC.)

- 7) When turning off the lamp, adjust the light intensity to the min and then turn off the power.

5. Photobiological safety of lamp and lamp systems.

CAUTION

This product is a light source which is classified as Risk Group 2 for blue light hazard and blue light small source hazard.

Possibly hazardous optical radiation emitted from this product.
Do not stare at operating lamp. May be harmful to the eyes.


< JCR15V150W5H/5 >

Risk Group 2 (Blue light hazard, Blue light small source hazard)
Other hazards: Exempt Group

IEC62471 explanatory

Caution : hot surface label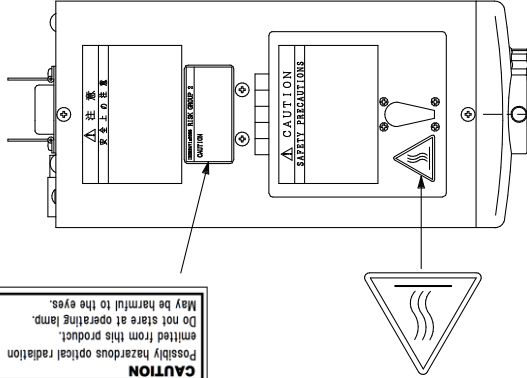

CAUTION

Because the cover stopper is hot, please be careful when you touch.

WARNING

Don't remove the heat absorbing filter and cold filter in the light source.
Removing the filters, the product is classified risk group 3.

When removing a heat absorbing filter and cold filter, the product is classified following risk group, because of the evaluation measurement of the IEC62471 safety.

< JCR15V150W5H/5 without filters >

Risk Group 3 (Active UV hazard)
Risk Group 2 (Blue light hazard, Blue light small source hazard, Retinal thermal hazard)
Risk Group 1 (Near UV hazard, IR radiation, IR radiation, eye hazard)
Other hazards: Exempt Group

< JCR12V100W10H/AL without filters >



Risk Group 3 (Active UV hazard)
Risk Group 2 (Blue light hazard, Blue light small source hazard, Retinal thermal hazard, IR radiation, eye hazard)
Other hazards: Exempt Group

- ① Power switch
 - This switch is used to turn on / off the power.
- ② Lamp burnout warning LED (Red)
 - This LED is illuminated when the lamp burns out or is not installed.
- ③ Light quantity control knob
 - This knob is used when controlling brightness of the lamp manually. A light quantity is increased by turning the knob in the clockwise direction.
- ④ Light guide socket
 - This is a socket to introduce a light from the halogen lamp built in this light source to the light guide.
- ⑤ Lamp house opening / closing cover
 - When you replace the lamp, loosen the lamp replacement hatch fixing screw to carry out replacement.
- ⑥ Suction ports
 - These are suction ports to allow the open air into this light source . For better ventilation, keep the suction ports 50mm or more away from an obstacle such as wall. If these ports are blocked, an light source internal temperature will increase and lead to a trouble.

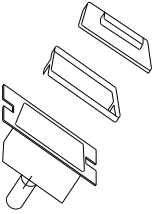
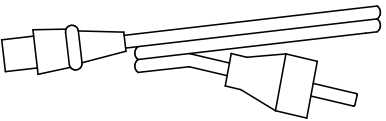
4. INSTALLATION

Do not use the product placed upside down or sideways down and the condition to cover ventilation ports. (suction, exhaust)
 Do not put the product on the position that a power supply-cord cannot be removed.
 Be sure to connect the protective conductor terminal of the power supply-cord to ground.
 The safety system of the product is off in the any system which the manufacturer do not authorize. In this case, the user has the responsibility of the safety on the system.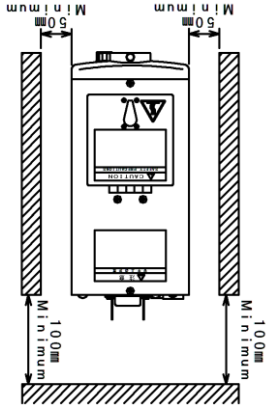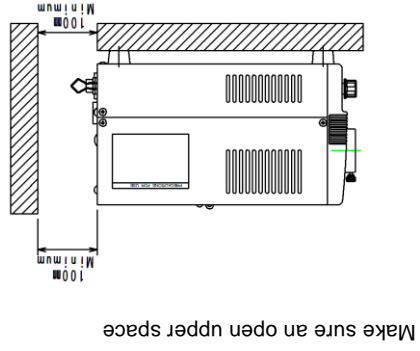

2. Items included the package


 <p>① Instruction manual</p>	 <p>② Plug for Remote Control (150CE, 150UE-B, 150UE-C)</p>	 <p>③ Power supply-cord (150UE-B: USA/Japan Type) (150CE, 150UE-C: European Type)</p>
--	--	--

CAUTION

Make sure to use the power supply-cord specified in "Specification." Using a wrong power supply-cord may result in malfunctions or fire. To prevent an electric shock, turn off the power switch (flip to OFF side) of the product whenever connecting or disconnecting the power supply-cord.

3. COMPONENT PARTS

3 - 1 Names of Component Parts

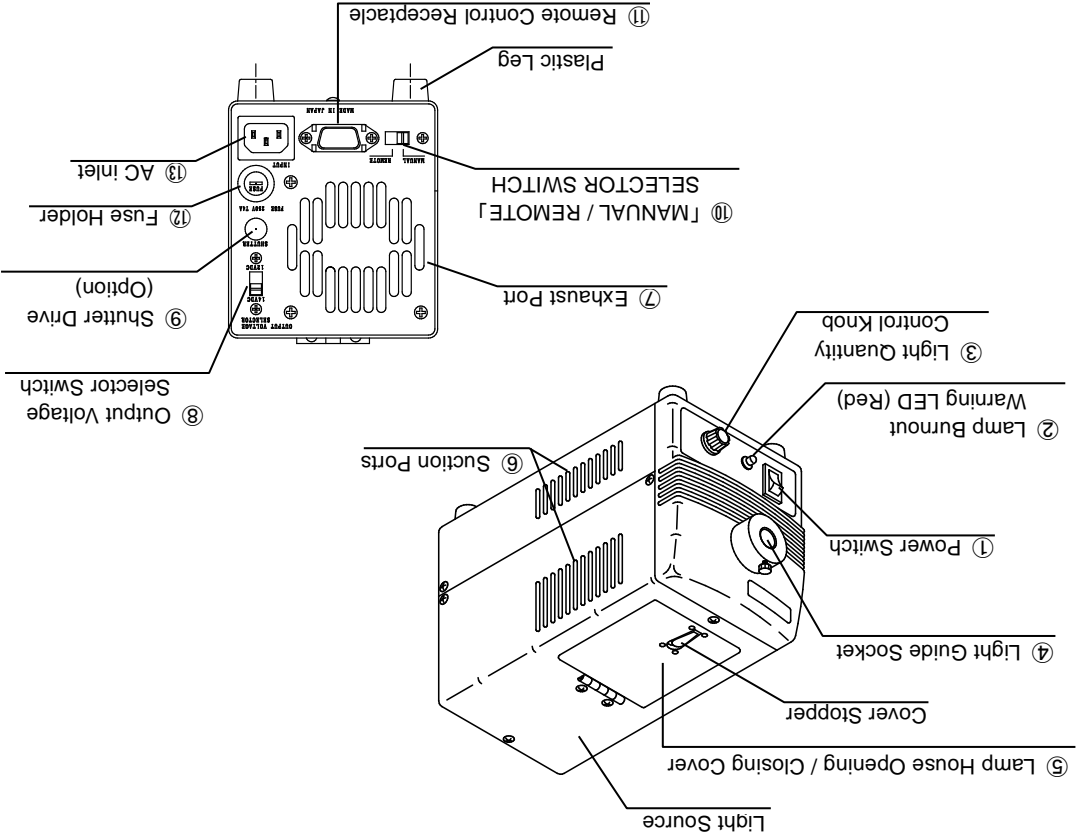LIGHT SOURCE

LUMINAR ACE

Model LA-150CE

LA-150E-B

LA-150E-C

INSTRUCTION MANUAL

Thank you for purchasing our product.

Prior to using it, please read the instruction manual thoroughly for safe and correct operation. Be sure to keep the instruction manual carefully at hand.

This light source is used for the illumination of image processing, microscope, science observation. This light source can not be installed in any system, and used with environment and installation which exceed the specification of product.

1. SAFETY PRECAUTIONS

PRECAUTIONS FOR SAFETY

This means that injury to the operator and damage to the equipment may result if these instructions are ignored or the equipment handled incorrectly.

- Watch out for High Temperature
While the product is being illuminated or immediately after it is turned off, never touch the lamp of the product or its turned surroundings, because they have a high temperature. (You may be burnt.)

- Watch out for Ignition
Never bring a combustible substance close to the illuminated product or its lamp. (It may cause a fire.)

- Watch out for Electric Shock
When you attach / detach a lamp, replace a fuse, or wire the cables, be sure to disconnect a power plug from a plug socket. (You may get an electric shock.)

- Do not look straight into the Lamp
Do not look at they illuminated lamp or a light coming through a light guide closely. (It may hurt your eyes or damage your eyesight.)

PRECAUTIONS FOR USE

- Do not move, vibrate, or give a shock to the product while it is being illuminated. (The lamp may burnout.)

- Do not use an unspecified lamp or plug socket. (A trouble may result.)

- Pay attention how a fan is running. Do not block a suction port or exhaust port or exhaust port while it is running. (A trouble may result.)

- Do not disassemble or re-engineer the product. (A trouble may result.)

- When replacing the lamp or plug socket, be sure to follow the instructions mentioned in REPLACEMENT OF MAINTENANCE PARTS in the instruction manual.

- When you turn on the lamp or while it is being illuminated, be sure to close a lamp house opening / closing cover.

- Avoid using the product in a place exposed to dust, oil, or iron powder, because it may shorten the service life of the lamp or damage the product.

- Do not install the product near a strong magnetic field or strong electric noise source. Malfunction may result from this.